pHY Marker

Code No. 3404A **Size:** **20 μ g**
Conc.: **0.2 - 0.8 μ g/ μ l**

Supplied Reagents:
6X Loading Buffer **1 ml**

* Please refer to the concentration of each lot in Certificate of Analysis supplied with the product.

Form:
10 mM Tris-HCl, pH 8.0
1 mM EDTA

Storage: -20°C
(Once opened, 6X Loading Buffer should be stored at room temperature.)
* 2 years from date of receipt under proper storage conditions.

Description:
This Marker is prepared by mixing pHY300PLK plasmid digested by restriction enzyme *Hind* III and pHY300PLK dimer plasmid digested by restriction enzyme *Hae* III.

Fragment	Size (bp)	Ratio * (%)
A	4,870	7.7
B	2,016	6.4
C	1,360	25.8
D	1,107	17.5
E	926	14.6
F	658	12.5
G	489	9.2
H	267	5.1
I	80	1.5

* Weight percentage of each fragment in total DNA

Usage:
Used as a DNA molecular size marker in agarose gel electrophoresis. This product can be radioactively labeled by the kinase exchange reaction using MEGALABEL™ (Cat. #6070).*

* Not available in all geographic locations. Check for availability in your area.

6X Loading Buffer (Store at RT after used):

36% Glycerol
30 mM EDTA
0.05% Bromophenol Blue
0.035% Xylene Cyanol

Add 1/5 volume of 6X Loading Buffer with DNA solution to apply on agarose gel electrophoresis. In case precipitates generated during the storage at room temperature, dissolve in warm bath before use.

Application example:

pHY Marker	X μ l (200 - 500 ng)
6X Loading Buffer	1 μ l
TE Buffer	up to 6 μ l

Run agarose gel electrophoresis using ex. 1% Agarose L03 [TAKARA] (Cat. #5003).

Perform staining with EtBr, or SYBR® Green I Nucleic Acid Gel Stain (Cat. #5760A/5761A).*

* Not available in all geographic locations. Check for availability in your area.

Manufactured by Yakult Co., Ltd.

SYBR is a registered trademark of Life Technologies Corporation. MEGALABEL is a trademark of Takara Bio Inc.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com. Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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pHY Marker	X μ l (200 ~ 500 ng)
6 \times Loading Buffer	1 μ l
TE Buffer	up to 6 μ l

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