

SeaKem® Gold Agarose

The fastest agarose available for separation of megabase DNA.

Introduction

SeaKem Gold Agarose is a unique, patented, very high gel strength, low EEO (≤ 0.05) standard gelling temperature agarose. This Genetic Technology Grade™ Agarose product was developed for rapid resolution of DNA and PCR[†] products between 1 kb and 50 kb by conventional electrophoresis.

Analytical Specifications

| | |
|----------------------------|-----------------------------|
| Gelling temperature (1.5%) | 36°C \pm 1.5°C |
| Melting temperature (1.5%) | $\geq 90^\circ\text{C}$ |
| Gel strength (1%) | $\geq 1,800 \text{ g/cm}^2$ |
| Gel strength (1.5%) | $\geq 3,500 \text{ g/cm}^2$ |

Applications

- Preparative DNA and RNA electrophoresis
- Analysis of Long PCR[†] reactions
- Separation and further manipulation of DNA $\geq 1,000$ bp
- Analytical electrophoresis of DNA and RNA
- Blotting of DNA and RNA

Suggested Agarose Concentrations

| Size Range (Base Pairs) | Final Agarose Concentration (%) 1X TAE Buffer |
|----------------------------|--|
| 5,000-50,000 | 0.3 |
| 1,000-20,000 | 0.5 |
| 800-10,000 | 0.8 |
| 400-8,000 | 1.0 |

† TBE Buffer is not recommended for separation of DNA $\geq 12,000$ bp.

Dye Mobility Table

Migration of double-stranded DNA in relation to Bromophenol Blue (BPB) and Xylene Cyanol (XC) in SeaKem Gold Agarose Gels.

| 1X TAE Buffer | | % Agarose | 1X TBE Buffer | |
|---------------|-------|--------------|---------------|-------|
| XC | BPB | | XC | BPB |
| 24,800 | 3,550 | 0.30 | 19,000 | 2,550 |
| 12,200 | 2,050 | 0.50 | 9,200 | 1,500 |
| 9,200 | 1,050 | 0.75 | 7,100 | 800 |
| 6,100 | 760 | 1.00 | 4,000 | 500 |
| 4,100 | 600 | 1.25 | 2,550 | 350 |
| 2,600 | 400 | 1.50 | 1,900 | 250 |
| 2,000 | 330 | 1.75 | 1,400 | 180 |
| 1,500 | 250 | 2.00 | 1,000 | 100 |

Precautions

Always wear eye protection when dissolving agarose and guard yourself and others against scalding solutions. Refer to Material Safety Data Sheet for additional safety and handling precautions.

Microwave Instructions for Agarose Preparation

1. Choose a beaker that is 2-4 times the volume of the solution.
2. Add room temperature 1X or 0.5X electrophoresis buffer and a stir bar to the beaker.
3. Slowly sprinkle in the agarose powder while the solution is rapidly stirred.
4. **Remove the stir bar if not Teflon® coated.**
5. Weigh the beaker and solution before heating.
6. Cover the beaker with plastic wrap.
7. Pierce a small hole in the plastic wrap for ventilation.
8. Heat the beaker in the microwave oven on **High** power until bubbles appear.
9. Remove the beaker from the microwave oven.
Caution: Any microwaved solution may become superheated and foam over when agitated.
10. **GENTLY** swirl the beaker to resuspend any settled powder and gel pieces.
11. Reheat the beaker on **HIGH** power until the solution comes to a boil.
12. **Hold at boiling point for 1 minute** or until all of the particles are dissolved.
13. Remove the beaker from the microwave oven.
14. **GENTLY** swirl the beaker to thoroughly mix the agarose solution.
15. After dissolution, add sufficient hot distilled water to obtain the initial weight.
16. Mix thoroughly.
17. Cool the solution to 50°C-50°C prior to casting.

Hot Plate Instructions for Agarose Preparation

1. Choose a beaker that is 2-4 times the volume of the solution.
2. Add room temperature electrophoresis buffer and a stir bar to the beaker.
3. Slowly sprinkle the agarose powder while the solution is rapidly stirred.
4. Weigh the beaker and solution before heating.
5. Cover the beaker with plastic wrap.
6. Pierce a small hole in the plastic wrap for ventilation.
7. Bring the solution to a boil while stirring.
8. Maintain gentle boiling until all the agarose is dissolved (approximately 10 minutes).
9. Add sufficient hot distilled water to obtain the initial weight.
10. Mix thoroughly.
11. Cool the solution to 50°C-60°C prior to casting.

Cambrex Bio Science Rockland, Inc.

www.cambrex.com/bioproducts

biotechserv@cambrex.com

Orders: 800-638-8174 Technical Service: 800-521-0390

Copyright © 2003 Cambrex Bio Science Rockland, Inc.



Ordering Information

| Catalog No. | Size |
|-------------|-------|
| 50150 | 125 g |
| 50152 | 25 g |

Related Products

AccuGENE® 10X TAE Buffer
GelStar® Nucleic Acid Gel Stain
Megabase DNA Standards
InCert® Agarose

†The PCR process may be covered by one or more third-party patents.

For *In Vitro* Diagnostic and Research Use.

For more information contact Technical Service at
(800) 521-0390 or visit our website at
www.cambrex.com

Cambrex Bio Science Rockland, Inc.

191 Thomaston Street
Rockland Maine 04841

To Order:

Phone: (800) 638-8174
Fax: (800)362-5552
E-mail: biocustserv@cambrex.com

Outside USA:

Phone: (207) 594-3400
Fax: (207) 594-3491

Technical Service:

Phone: (800) 521-0390
Fax: (800) 362-1133
E-mail: biotechserv@cambrex.com
Homepage: www.cambrex.com

Cambrex Bio Science Copenhagen, ApS.

Risingevej 1
DK-2665 Vallensbaek Strand
Denmark
Phone: 45 43 56 74 00
Fax: 45 43 56 74 03

Cambrex Bio Science Wokingham, Ltd.

1 Ashville Way
Wokingham, Berkshire RG41 2PL
United Kingdom
Phone: 44 118 979 5234
Fax : 44 118 979 5231

AccuGENE, GelStar, Genetic Technology Grade and InCert Agarose are trademarks of CBM Intellectual Properties inc., SeaKem and GelStar are trademarks of FMC Corporation. Teflon is a trademark of E.I. duPont de Nemours and Co.

© 2004 Cambrex Bio Science Rockland, Inc.
All rights reserved.

18106-0604-6

Cambrex Bio Science Rockland, Inc.
www.cambrex.com/bioproducts
biotechserv@cambrex.com

Orders: 800-638-8174 Technical Service: 800-521-0390
Copyright © 2003 Cambrex Bio Science Rockland, Inc.

