10XPCR Buffer, MgCl<sub>2</sub>, dNTP Mixture for PCR

# TaKaRa Ex Taq™

Code No. RR01AM

Size: 250 units

Shipping at -20°C

Stored at -20°C

Contents : TaKaRa Ex Taq™ (TaKaRa Cat.# RR001)  
10X Ex Taq™ Buffer (Mg<sup>2+</sup> free)  
25 mM MgCl<sub>2</sub>  
dNTP Mixture

Lot No.

Conc. : units/μl

Volume : μl

Expiry Date :

#### Storage Buffer:

20 mM	Tris-HCl (pH8.0)
100 mM	KCl
0.1 mM	EDTA
1 mM	DTT
0.5%	Tween® 20
0.5%	Nonidet P-40®
50%	Glycerol

**Unit definition:** One unit is the amount of the enzyme that will incorporate 10 nmol of dNTP into acid-insoluble products in 30 minutes at 74°C with activated salmon sperm DNA as the template-primer.

#### Reaction mixture for unit definition:

25 mM	TAPS (pH 9.3 at 25°C)
50 mM	KCl
2 mM	MgCl <sub>2</sub>
1 mM	2-mercaptoethanol
200 μM	each dATP, dGTP, dTTP
100 μM	[α- <sup>32</sup> P]-dCTP
0.25 mg/ml	activated salmon sperm DNA

**Purity:** Nicking activity, endonuclease and exonuclease activity were not detected after the incubation of 0.6 μg of supercoiled pBR322 DNA, 0.6 μg of λ DNA or 0.6 μg of λ -Hind III digest with 10 units of this enzyme for 1 hour at 74°C.

#### Applications:

For DNA amplification by Polymerase Chain Reaction (PCR).

**PCR products :** As most PCR products amplified with TaKaRa Ex Taq™ have one A added at 3'-termini, the obtained PCR product can be directly used for cloning into T-vector. Also it is possible to clone the product in blunt-end vectors after blunting and phosphorylation of the end.

**PCR test :** Good performance of DNA amplification by Polymerase Chain Reaction (PCR) was confirmed by using λ DNA as the template (amplified fragment : 20 kbp).

Good performance of DNA amplification of β-globin gene by PCR was also confirmed by using human genome DNA as the template (amplified fragment : approx 3 kbp).

#### General reaction mixture for PCR (total 50 μl)

TaKaRa Ex Taq™ (5 units/μl)	0.25 μl
10X Ex Taq™ Buffer (Mg <sup>2+</sup> free)	5 μl
MgCl <sub>2</sub> (25 mM)	4 μl
dNTP Mixture (2.5 mM each)	4 μl
Template	< 500 ng
Primer 1	0.2 ~ 1.0 μM
Primer 2	0.2 ~ 1.0 μM
Sterilized distilled water	up to 50 μl

#### Supplied 10X Ex Taq™ Buffer (Mg<sup>2+</sup> free)

Supplied Size	: 1 ml/vial
Storage	: -20°C

#### Supplied 25 mM MgCl<sub>2</sub>

Supplied Size	: 1 ml/vial
Storage	: -20°C

#### Supplied dNTP Mixture

Mixture of dNTP, ready for use in Polymerase Chain Reaction (PCR) without dilution.

Supplied Size	: 800 μl/vial
Concentration	: 2.5 mM of each dNTP
pH	: pH7~ 9
Form	: Solved in water (sodium salts)
Purity	: ≥ 98% for each dNTP
Storage	: -20°C

#### < Cool Start Method >

"Cool Start Method", enables to minimize the amplification of non-specific band in PCR and achieves more accurate amplification. This is a simpler method without need for special enzyme nor additional reagents.\* Higher reaction specificity can be achieved by combining Hot Start PCR techniques with Taq Antibody (Code.9002A) and Cool Start method.

#### Protocol of Cool Start Method

- 1) Keep all reagents on ice until use.
  - 2) Prepare the reaction mixture on ice.\*\*  
\*The adding order of reagents dose not influence on results.  
\*\*The result will not be affected even when the mixture is left on ice 30 min. before thermal cycling.
  - 3) Set a thermal cycler ready to start with the designated program.\*\*  
\*\*\*No need to change PCR conditions especially for Cool Start.
  - 4) Set the tubes in a thermal cycler and start thermal cycling immediately.
- \* JAPAN Patent 2576741 for Cool Start Method is owned by SHIMAZU CORPORATION.

#### Note

For research use only. Not for use in diagnostic or therapeutic procedures.

Produced by TAKARA BIOTECHNOLOGY (DALIAN) CO.,LTD.

U.S. Patent 5,436,149 for LA Technology is owned by TAKARA BIO INC.

#### NOTICE TO PURCHASER: LIMITED LICENSE

A license under the foreign counterparts of U.S. Patents Nos. 4,683,202, 4,683,195 and 4,965,188 owned by F.Hoffmann-La Roche Ltd. and U.S. Patent No. 5,075,216 and its foreign counterpart, owned by Roche Molecular Systems, Inc. and F.Hoffmann-La Roche Ltd. for use in research and development, has an up-front fee component and a running-royalty component. The purchase price of this product includes limited, nontransferable rights under the running-royalty component to use only this amount of the product to practice the polymerase chain reaction (PCR) and related processes described in said patents where such processes are covered by patents solely for the research and development activities of the purchaser when this product is used in conjunction with a thermal cycler whose use is covered by the up-front fee component. Rights to the up-front fee component must be obtained by the end user in order to have a complete license to use this product in the PCR process where the process is covered by patents. These rights under the up-front fee component may be purchased from Applied Biosystems or obtained by purchasing an authorized thermal cycler. No right to perform or offer commercial services of any kind using PCR, where the process is covered by patents, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is hereby granted by implication or estoppel. Further information on purchasing licenses to practice the PCR process where the process is covered by patents may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94044 or the Licensing Department, Roche Molecular Systems, Inc., 1145 Atlantic Avenue, Alameda, California 94501.

10 × PCR Buffer, MgCl<sub>2</sub>, dNTP Mixture for PCR

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