Code RR260A

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Takara

***Kudoa septempunctata* (18S rRNA)**

qPCR Detection Kit

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v201302Da

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NOTICE TO PURCHASER: LIMITED LICENSE

[A6] PCR Notice

Use of this product is covered by one or more of the following US patents and corresponding patent claims outside the us: 5,804,375,5,538,848,5,723,591,5,876,930,6,030,787 and 6,258,569. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product solely in Environmental Testing, Food Testing, Industrial Microbiology, including reporting results of purchaser's activities for a fee or other commercial consideration, and also for the purchaser's own internal research. No right under any other patent claim is conveyed expressly, by implication, or by estoppel. Further information on purchasing licenses may be obtained from the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

[L15] Hot Start PCR

Licensed under U.S. Patent No. 5,338,671 and 5,587,287 and corresponding patents in other countries.

[L53a] Rox Reference Dye (Application Field)

Use of this product is covered by one or more of the following US patents and corresponding patent claims outside the US: 5,928,907. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product solely in Environmental Testing, Industrial Microbiology, including reporting results of purchaser ' s activities for a fee or other commercial consideration, and also for the purchaser's own internal research. No right under any other patent claim is conveyed expressly, by implication, or by estoppel. Further information on purchasing licenses may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

[M40] Thermostable RNaseH

This product is covered by the claims of U.S. Patent No. 7,422,888 and its foreign counterpart patent claims.

[M57] LA Technology

This product is covered by the claims 6-16 of U.S. Patent No. 5,436,149 and its foreign counterpart patent claims.

[M82] Tli RNaseH Plus

This product is the subject of the pending JP patent application.