

Stellar™ Competent Cells Protocol-at-a-Glance

(PT5055-2)

Stellar Competent Cells are an *E.coli* HST08 strain that provides high transformation efficiency paired with blue-white screening capability when used with pUC plasmid vectors.

Genotype

F⁻, endA1, supE44, thi-1, recA1, relA1, gyrA96, phoA, Φ80d lacZΔ M15, Δ (lacZYA - argF) U169, Δ (mrr - hsdRMS - mcrBC), ΔmcrA, λ-

Please Read Before Proceeding with Transformation

1. Use no more competent cells than necessary. Transport cells on dry ice/ethanol.
2. You may use 1.5-ml microcentrifuge tubes instead of 14-ml round-bottom tubes for transformation, but this may reduce efficiency.
3. When transforming 100 μl of competent cells, do not use more than 10 ng of purified sample DNA. If you use more than 10 ng of DNA, transformation efficiency may decrease.
4. If you change the amount of competent cells, or types of tubes used, it might be necessary to reevaluate the optimal conditions. For example, when using 1.5-ml microcentrifuge tubes, heat shock for 60 seconds at 42°C (see Transformation Protocol Step 5, below).
5. When adding X-Gal to medium, do so as follows:
Add 20 mg/ml X-Gal (dissolved in dimethylformamide) into 200 μl/100 ml agar medium.
6. Do not refreeze competent cells once thawed. If necessary, freeze the cells in dry ice and stock at -70°C. However, the transformation efficiency may decrease more than one order of magnitude.

Transformation Protocol

1. Thaw Stellar Competent Cells in an ice bath just before use.
2. After thawing, mix gently to ensure even distribution, then move 100 μl of competent cells into a 14-ml round-bottom tube (falcon tube). Do not vortex.
3. Add no more than 10 ng of DNA for transformation.
4. Place tubes on ice for 30 min.
5. Heat shock the cells for exactly 45 sec at 42°C.
6. Place tubes on ice for 1–2 min.
7. Add SOC medium to bring the final volume to 1 ml. SOC medium should be warmed to 37°C before using.
8. Incubate by shaking (160–225 rpm) for 1 hr at 37°C.
9. Plate an appropriate amount of culture on selective medium.*
***Note:** For a plate with a diameter of 9 cm, plate 100 μl. Plating is accomplished by spreading cells on selective medium [e.g., LB agar + Ampicillin(50–100 μg/ml)]. The medium should also contain X-gal (40 μg/ml) for plasmids that permit blue/white screening of transformants.
10. Incubate overnight at 37°C.



Clontech

United States/Canada
800.662.2566

Asia Pacific
+1.650.919.7300

Europe
+33.(0)1.3904.6880

Japan
+81.(0)77.543.6116

Clontech Laboratories, Inc.
A Takara Bio Company
1290 Terra Bella Ave.
Mountain View, CA 94043
Technical Support (US)
E-mail: tech@clontech.com
www.clontech.com

Notice to Purchaser

Clontech products are to be used for research purposes only. They may not be used for any other purpose, including, but not limited to, use in drugs, *in vitro* diagnostic purposes, therapeutics, or in humans. Clontech products may not be transferred to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service to third parties without written approval of Clontech Laboratories, Inc.

Clontech, the Clontech logo and all other trademarks are the property of Clontech Laboratories, Inc., unless noted otherwise. Clontech is a Takara Bio Company. ©2009 Clontech Laboratories, Inc.

(PR983324; published August 2009)