pUC 18 DNA

Code No. 3218

Size: 25 μ g

Conc.: 0.5 μ g/ μ l

* 2 years from date of receipt under proper storage conditions.

Description:

This DNA is a plasmid vector suitable for DNA sequencing by dideoxy method. This DNA may allow successful cloning of DNA fragment longer than the case of M13 phage vector used. Since it has multi-cloning sites on *lacZ'* region, any DNA insertion into this vector can be easily verified using plates containing IPTG and X-Gal. Moreover, expression of the cloned DNA can be performed by using of *lac* promoter. M13 primers are available for DNA sequencing. Using Kilo-Sequence Deletion Kit (Cat. #6030), DNA sequencing of the cloned DNA with kilo-bases can also be performed.


Form: 10 mM Tris-HCl, pH 8.0
1 mM EDTA

Storage: -20°C

Preparation: Purification of cccDNA by CsCl-EtBr ultracentrifugation

Chain length: 2,686 bp

Cloning site of pUC18 :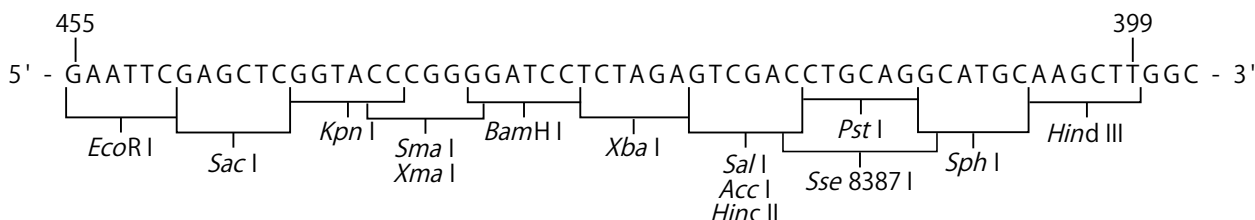

Quality Control Data:

Please see the Certificate of Analysis (CoA) for each lot. You can download the CoA on Takara Bio website.

Usage:


- Cloning of a target gene and its expression using *lac* promoter.
- DNA sequencing by using M13 primers.
- Determination of long DNA sequences using Deletion Kit for Kilo-Sequencing (Cat. #6030).

Reference:

- 1) Yanisch-Perron C, Vieira J, and Messing J. *Gene*. (1985) **33**: 103-119.
- 2) Messing J. *Methods in Enzymology*. (1983) **101**: 20-78.

GenBank:

Enty Name : SYNPU18CV
Accession No : L09136

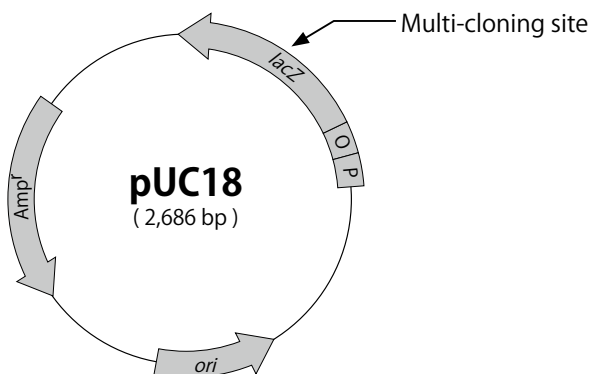Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com. Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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