CLEARLY Stained Protein Ladder

Code No. 3454A Size: 500 μ l

Description :

CLEARLY Stained Protein Ladder is a mixture of 11 pre-stained recombinant highly purified proteins. When this ladder is analyzed by SDS-PAGE, it is resolved into clearly identifiable visible bands from 10 to 210 kDa. The 30 and 70 kDa bands have quadruple the mass of the other bands and serve as reference indicators.

This ladder is pre-mixed with loading buffer and ready to load. Pre-heating is not required. 10 μ l of this ladder is loaded per lane of SDS-PAGE minigel for standard use. In this case, this ladder is sufficient for 50 lanes.


This ladder is useful for monitoring of protein migration during SDS-PAGE and verifying transfer efficiency at western blotting.*

- * Because the covalent coupling of the dye to the proteins affects their electrophoretic behavior on SDS-PAGE gels relative to unstained proteins, this ladder is not recommended for use as a precise size standard when performing SDS-PAGE to calculate the molecular weight of a protein of interest.

Storage: -20°C

- * This ladder is stable for 6 months if stored at 4°C.
- * This ladder is stable after 50 times of freeze-thaw cycles.

< Electrophoresis Result (12% SDS-PAGE) >

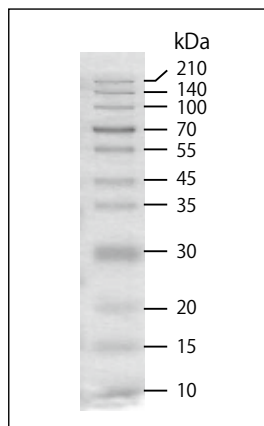No staining

Application Example :

1. Thaw CLEARLY Stained Protein Ladder at room temperature. Confirm the ladder is completely dissolved. Pre-heating is not required.
2. Load 10 μ l per lane of SDS-PAGE minigel. Run SDS-polyacrylamide gel electrophoresis.

Note :

It is recommended to use 5 - 20% gradient SDS-polyacrylamide gel for electrophoresis. If the concentration of gel is too low, the mobility of bands of low molecular weight is faster than the mobility of BPB. If the concentration of gel is too high, high molecular weight bands may not separate well or they may not migrate.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc.

If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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