

Cat. # 9021

For Research Use

TAKARA

***E. coli* HB101 Electro-Cells**

Product Manual

v202010Da

Table of Contents

I.	Description.....	3
II.	Components	3
III.	Storage	3
IV.	Protocol.....	4
V.	Quality.....	5
VI.	Genotype.....	5
VII.	Cell Density.....	5
VIII.	References.....	5
IX.	Related Products	5

I. Description

E. coli HB101 are specially prepared by Takara Bio to be best appropriate for electroporation method. Electroporation method is used to transfer DNA into a cell by breaking cytoplasmic membrane by high voltage pulse. As this electro-cells offer high transformation efficiency and good reproducibility, it is especially useful in transferring small amount of sample into an *E. coli* in less time.

E. coli HB101 Electro-Cells can be used for preparation of DNA library or subcloning of recombinant plasmid, that does not require blue/white selection.

II. Components

<i>E. coli</i> HB101 Electro-Cells	50 μ l x 10
pBR322 plasmid (10 pg/ μ l)	10 μ l
SOC Medium*	1 ml x 10

* SOC Medium	2%	Tryptone
	0.5%	Yeast extract
	10 mM	NaCl
	2.5 mM	KCl
	10 mM	MgSO ₄
	10 mM	MgCl ₂
	20 mM	Glucose

III. Storage

-80°C

Note : Store at -80°C or lower. If the storage temperature is not maintained consistently, the transformation efficiency will be reduced. You may determine the transformation efficiency of stored cells by using the included pBR322 control.
Do not store in liquid nitrogen.

IV. Protocol

- (1) Thaw 50 μ l of *E. coli* HB101 Electro-Cells in an ice bath just before use.
- (2) Add 1 - 2 μ l of DNA solution* into the thawed cell suspension.
 - * When sample DNA solution contains salt, dilute with TE buffer or sterile purified water. Or desalting by ethanol precipitation is recommended (≤ 10 ng is recommended).
- (3) Transfer the mixture of cells and DNA to a cold 0.1 cm electroporation cuvette.
- (4) After applying pulse*, immediately add 1 ml of SOC Medium (precooled in an ice bath).
 - * Takara Bio uses BIO-RAD MicroPulser and the electrical condition is 1.8 kV. In the case of BIO-RAD Gene Pulser, standard electrical conditions are 200 Ω , 25 μ F, 1.8 kV.
- (5) Transfer into a 14 ml round-bottom tube (CORNING Code: 352059 or 352057, etc.). Incubate by shaking (160 - 250 rpm) for 1 hour at 37°C.
- (6) Plate on selective media.*
 - * Plate no more than 100 μ l for a ϕ 9 cm plate. If necessary, dilute the culture with the same medium as used in step (4).
- (7) Incubate overnight at 37°C.

【 Read these before use 】

1. Place a vial of electro-cells in a dry ice / EtOH bath immediately upon removal from -80°C freezer. Keep cells in bath until you are ready to proceed.
2. When using 50 μ l of electro-cells, apply high-purified sample DNA in less than 10 ng. If not, transformation efficiency might decrease.
3. When changing an experiment scale, optimum condition should be considered.
4. When transferring high molecular weight DNA (> 7 kb), transformation efficiency may decrease.
5. Use TE buffer for sample DNA preparation. High salt concentration in sample DNA solution may decrease transformation efficiency.
6. L-broth or ϕ b-broth can be used instead of SOC Medium. In this case, lower efficiency might be obtained.

<u>L-broth</u> :	<u>Ingredient</u>	<u>per liter water</u>
	Tryptone	10 g
	Yeast extract	5 g
	NaCl	5 g

Adjust to around pH 7.5 with 1 N NaOH and autoclave.

<u>ϕ b-broth</u> :	<u>Ingredient</u>	<u>per liter water</u>
	Tryptone	20 g
	Yeast extract	5 g
	MgSO ₄ ·7H ₂ O	5 g

Adjust to around pH 7.5 with 1 N KOH and autoclave.

<u>L-plates</u> :	<u>Ingredient</u>	<u>per liter water</u>
	Tryptone	10 g
	Yeast extract	5 g
	NaCl	5 g

Adjust to around pH 7.5 with 1N NaOH, add agar to be 1.5% and autoclave.

7. Once the Electro-Cells have been thawed, refreezing for storage is not recommended. If this is unavoidable, flash freeze the cells on dry ice/ethanol and store them promptly at -80°C. However, the transformation efficiency will be lowered by at least one order of magnitude.

V. Quality

10 pg of pBR322 was transformed and selected by plating to Amp⁺ selective media. Transformation efficiency : > 5 x 10⁸ cfu/ μg pBR322

VI. Genotype

E. coli HB101: *supE44*, Δ (*mcrC-mrr*), *recA13*, *ara-14*, *proA2*, *lacY1*, *galK2*, *rpsL20*, *xyl-5*, *mtl-1*, *leuB6*, *thi-1*.

VII. Cell density

>1 x 10¹⁰ bacteria/ml

VIII. References

- 1) Dower W J, Miller J F, and Ragsdale C W. *Nucl Acids Res.* (1988) **16**: 6127.
- 2) Bottger E C. *Biotechniques.* (1988) **6**: 878.

IX. Related products

E. coli HB101 Competent cells (Cat. #9051)
pBR 322 DNA (Cat. #3050)

NOTE: This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc.

Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc.

If you require licenses for other use, please contact us by phone at +81 77 565 6972 or from our website at www.takara-bio.com.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.
