6X Loading Buffer

Code No. 9156 **Size:** **1 ml x 10**

* 2 years from date of receipt under proper storage conditions.

Description:

This product is Loading Buffer for agarose gel electrophoresis. It is the same composition as 6X Loading Buffer supplied with DNA Digested Markers (λ -*Hind* III digest, etc) and DNA Ladder Markers (100 bp DNA Ladder, etc).

Composition :

36%	Glycerol
30 mM	EDTA
0.05%	Bromophenol Blue
0.035%	Xylene Cyanol

Storage: Room temperature

Applications:

For agarose gel electrophoresis.

Instruction:

Add 1/5 volume of 6X Loading Buffer with DNA solution to apply on agarose gel electrophoresis. In case precipitates generated during the storage at room temperature, dissolve in warm bath before use.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc.

If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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